
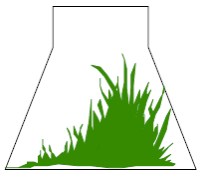


LILY MULTIPLICATION KIT

Product No. L577

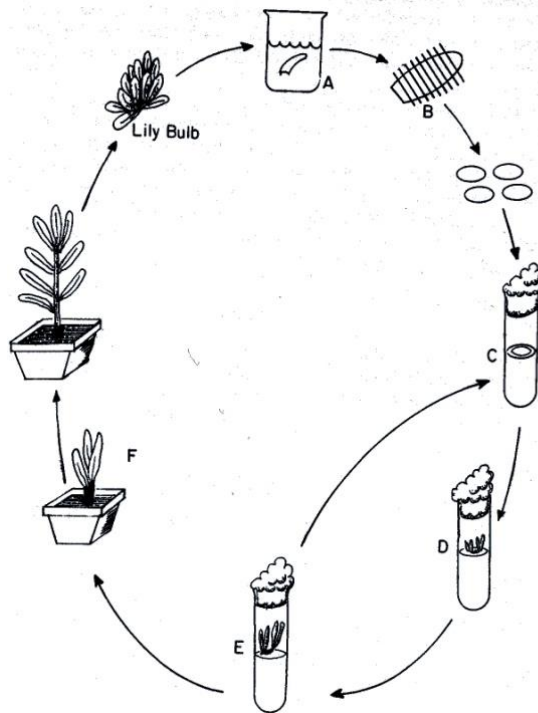

PhytoTechnology Laboratories®

Mailing Address: P.O. Box 12205, Shawnee Mission, KS 66282-2205

Phone: 1-888-749-8682 (1-913-341-5343 *Outside the USA & Canada*)

Fax: 1-888-449-8682 (1-913-341-5442 *Outside the USA & Canada*)

Visit our Web Site at <http://www.phytotechlab.com>

©2013 PhytoTechnology Laboratories®



TABLE OF CONTENTS

KIT COMPONENTS.....	2
MATERIALS REQUIRED BUT NOT PROVIDED.....	3
INTRODUCTION.....	3
MEDIA PREPARATION	3
STERILIZATION OF MEDIA	4
MEDIA STERILIZATION TIMES	4
CULTURE PROCEDURES.....	5
APPROXIMATE SCHEDULE.....	5
LITERATURE CITED	5
MEDIA FORMULATIONS.....	6
STOCK SOLUTION AND MEDIA PREPARATION LOG	7
NOTES.....	8

KIT COMPONENTS

Product No.	Product Description	1 Each
	Box	1
	Instruction Manual	1
C215 – 10 ea	Culture Containers	1
F951 – 1 ea	Forceps, 8”	1
S963 – 1 ea	Scalpel Handle, No.3	1
S971	Scalpel Blades, No. 11	2
P334 – 1 roll	pH Strips, 4.5 – 7.5	1
D940 – 20 ea	Petri Dishes	1
V886 – 15 mL	Vinegar	1
S803 – 25 g	Sodium Bicarbonate (Baking Soda)	1
P068	Pipette, Plastic Transfer	2
M401 – 1L	Murashige & Skoog Modified Basal Medium (w/ BA)	4
M508 – 1L	Murashige Fern Multiplication Medium	4
M555 – 1L	Murashige & Skoog Modified Multiplication Medium (w/ Kinetin)	4
S391 – 500 g	Sucrose	1
A296 – 9 g	Agar	12
Optional at an additional charge	Lily Bulb (NOT included in kits sent outside the continental USA)	1

MATERIALS REQUIRED BUT NOT PROVIDED

1. Beakers/containers: two 250-mL and one 500 mL
2. Media preparation container
3. 10% chlorine bleach solution supplemented with a few drops of Tween-20 (Product No. 720)
4. 1000 mL of sterile distilled water (Product No. W783)
5. Tissue culture water (distilled or deionized)
6. 70% Isopropyl alcohol
7. Bunsen or alcohol burner (Product No. B966 or B876, respectively)

INTRODUCTION

Hussey studied the effect of the cytokinin 6-benzylaminopurine (BA) on the release of axillary buds of *Lilium pyrenacium* Govan and an Asiatic hybrid using adventitious buds from stems as the explant. Optimal production of buds (i.e., 1 to 5 laterals) was promoted by BA concentrations of 2.0-8.0 mg/liter on a Murashige and Skoog (MS) medium. Hussey later found that shoot proliferation in *L. longiflorum* and *L. pyrenacium* was promoted by the use of a cytokinin plus an auxin when pieces of leaves were used as the explant source. Stem sections cultured on a MS medium supplemented with 2.0 mg/liter Indole-3-acetic acid (IAA) and 0.5 mg/liter BA also produced approximately 10 bulblets per section.

Steinberg et al. obtained optimal bulblet formation from leaf blades of *L. longiflorum* 'Nellie White' when cultured on a MS medium supplemented with 10.0 mg/liter Naphthaleneacetic acid (NAA) and 1.0 mg/liter kinetin. This combination, however, was the least effective in promoting shoot formation. Initiation of one to three bulbs occurred at the base of each leaf.

The purpose of this kit is to demonstrate vegetative propagation and the effects of the position that an explant is taken from bulb and the role it may play in shoot multiplication. This kit will also examine the effects of vitamins on shoot proliferation.

MEDIA PREPARATION

Powdered media are extremely hygroscopic and must be protected from atmospheric moisture. If possible, the entire contents of each package should be used immediately after opening. Media stored at 2 to 6°C and tightly sealed should last 2-3 years. Preparing the medium in a concentrated form is not recommended as some salts in the medium may precipitate. The basic steps for preparing the culture medium are listed below:

1. Measure out approximately 90% of the desired final volume of tissue culture grade water, e.g. 900 mL for a final volume of 1000 mL. Select a container twice the size of the final volume.
2. While stirring the water, add the powdered medium and stir until completely dissolved.
3. Rinse the container that the medium was packaged in with a small volume of tissue culture grade water to remove traces of the powder. Add to the solution in Step 2.
4. Add agar while stirring; it will not dissolve but should disperse into a uniform suspension.
5. Add 6-9 g/L of agar to all media. Add 30 g/L of sucrose to M508; it is already contained in M401 and M555. Aside from these supplements, the media provided in this kit are complete and typically do not require other supplements.
6. Add additional tissue culture grade water to bring the medium to the final volume.

7. While stirring, determine the pH using the pH Strips (Product No. P334). If necessary, adjust the medium to the desired pH using the baking soda to raise the pH or vinegar to lower the pH. A pH of 5.6 to 5.8 is typically recommended for most plants, including fern. Alternatively, the pH can be adjusted by using dilute potassium hydroxide or sodium hydroxide solution to raise the pH and dilute hydrochloric (muriatic) acid to lower the pH of the medium.
8. While stirring, heat the solution to nearly boiling to melt the agar in the medium.
9. Dispense the medium into the culture vessels before or after autoclaving as indicated below:
 - a. The Petri dishes (Product No. D940) included in this kit are sterile and cannot be autoclaved. They will melt if heated in an autoclave (or pressure cooker). Medium to be dispensed in Petri dishes must be sterilized and partially cooled before pouring it in the dishes.
 - b. The culture vessels (Product No. C215) are autoclavable. Media should be dispensed in these vessels prior to sterilization in an autoclave or pressure cooker. The lids of these culture vessels C215 should not be tightly sealed during sterilization to allow for proper steam and pressure penetration.
10. Sterilize the medium in a validated autoclave or pressure cooker at 1 kg/cm², 121 °C (15 psi, 250 ° F), for the time period described under “Sterilization of Media” below.
11. Allow medium to cool prior to use.

STERILIZATION OF MEDIA

Plant tissue culture media are generally sterilized by autoclaving at 121 °C and 1.05 kg/cm² (15 psi). This high temperature not only kills bacteria and fungi, but also their heat-resistant spores. Media can be sterilized in either an autoclave or pressure cooker with similar results. The time required for sterilization depends upon the volume of medium in the vessel. The minimum times required for sterilization of different media volumes are listed below. It is advisable to dispense medium in small aliquots whenever possible as many media components are broken down by prolonged exposure to heat.

MEDIA STERILIZATION TIMES

Volume of Medium per Vessel (mL)	Minimum Autoclaving Time (min.)
25	15-20
50	25
100	28
250	31
1000	40
2000	48
4000	63

Please Note: Minimum Autoclaving Time includes the time required for the liquid volume to reach the sterilizing temperature (121 °C) and 15 minutes at 121 °C (Burger, 1988). Times may vary due to differences in autoclaves. Validation with your autoclave or pressure cooker is recommended.

CULTURE PROCEDURES

1. Wipe down all surfaces of the transfer hood or work area with 70% isopropyl alcohol. Allow the hood to run for 15 min before beginning transfer operations. Place all the materials listed in the previous section under the hood. Place scalpels and forceps in a 250-ml beaker containing about 100 ml of 70% isopropyl alcohol.
2. Remove outer layer of damaged scales from the lily bulbs and discard. Under running water, begin removing the rest of the scales making certain to remove all soil; then place the scales in a 250ml beaker containing tap water. Remove the scales down to the point where they become quite narrow and small, then discard the rest of the bulb. Place the nylon mesh over the beaker and place under running tap water so that the water gently agitates the scales. Adding a small amount of detergent may be helpful in the cleaning process. Allow the scales to rinse in this manner for 5-10 min. Once the scales have been washed thoroughly, pour off all water, remove the nylon mesh, and transfer the beaker of scales to the laminar flow hood.
3. Pour enough 95% ethanol over the scales to cover them, swirl for 3-4 min, and decant the ethanol. Next pour the bleach solution over the scales and sterilize for 15 min before decanting off the solution. Rinse the scales three times in sterile distilled water with each rinse lasting approximately 1 min.
4. Place the culture vessels containing the media in the hood/work area.
5. All tools which now contact the tissue should be sterilized in alcohol and then flamed to remove any alcohol.
6. Transfer several sterilized scales to a sterile Petri dish. Section the scales into pieces about 4 mm wide. Note whether each section comes from the base or tip of the scale. After sectioning, transfer several sections to each culture vessel. Note whether each section comes from the base or tip of the scale. Place sections from both the tip and base on the medium. Once all cultures have been completed, place them in low light (e.g., fluorescent light) at 25 °C.
7. Once shoots have developed they can be subcultured (individually transferred) onto fresh medium for continued multiplication or removed and planted in potting soil.

APPROXIMATE SCHEDULE

Event	Timing
Isolation of fresh explants	Day 0
First appearance of nodules (organogenesis)	Day 15
Noticeable bulblet formation	Day 30
First subculture	Day 60 - 90
Transfer to soil	Day 90+ (When plantlets are large enough to handle)

LITERATURE CITED

- Burger, D.W. 1988. Guidelines for autoclaving liquid media used in plant tissue culture. HortScience 23:1066-1068.

MEDIA FORMULATIONS

All components express in mg/L	Murashige & Skoog Modified Basal Medium (w/ BA)	Murashige Fern Multiplication Medium	Murashige & Skoog Modified Multiplication Medium w/Kinetin
COMPONENT	M401	M508	M555
Ammonium Nitrate	1650	1650	1650
Boric Acid	6.2	6.2	6.2
Calcium Chloride, Anhydrous	332.2	333	332.2
Cobalt Chloride·6H ₂ O	0.025	0.025	0.025
Cupric Sulfate·5H ₂ O	0.025	0.025	0.025
Na ₂ EDTA·2 H ₂ O	37.26	-	37.26
Ferric Sodium EDTA	-	36.7	-
Ferrous Sulfate·7H ₂ O	27.8	-	27.6
Magnesium Sulfate	180.7	181	180.7
Manganese Sulfate·H ₂ O	16.9	16.9	16.9
Molybdc Acid (Sodium Salt)·2H ₂ O	0.25	0.25	0.25
Potassium Iodide	0.83	0.83	0.83
Potassium Nitrate	1900	1900	1900
Potassium Phosphate, Monobasic	170	170	170
Sodium Phosphate Monobasic	-	255	148
Zinc Sulfate·7H ₂ O	8.6	8.6	8.6
Adenine Hemisulfate	-	-	80
6-Benzylaminopurine (BA)	1	-	-
Glycine (Free Base)	2	-	-
Kinetin	-	2	1
myo-Inositol	100	100	100
a-Naphthaleneacetic Acid	0.1	0.1	0.1
Nicotinic Acid (Free Acid)	0.5	-	-
Pyridoxine·HCl	0.5	-	-
Sucrose	30000	-	30000
Thiamine·HCl	0.4	0.4	0.4
Grams of powder to prepare 1 liter	34.44	4.66	34.66

STOCK SOLUTION AND MEDIA PREPARATION LOG

Product Number: _____ Medium: _____

Lot Number: _____ Prepared By/ Date: _____

Volume to Prepare: _____ Autoclave Sterilization Time: _____

pH Desired: _____ Actual Final pH: _____

Instructions: Complete the table with all components of the stock solution or medium to be prepared, including the product number, lot number, and grams/batch. As each component is weighed record the actual weight on the sheet. Check off each component after it is added to the solution/medium.

Component	Product Number	Lot Number	Grams/ Batch	Actual Weight	Added <input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>
					<input type="checkbox"/>

Instructions/ Comments:

Species/Tissue Cultured: _____

NOTES

PhytoTechnology Laboratories®

Mailing Address: P.O. Box 12205, Shawnee Mission, KS 66282-2205

Phone: 1-888-749-8682 (1-913-341-5343 *Outside the USA & Canada*)

Fax: 1-888-449-8682 (1-913-341-5442 *Outside the USA & Canada*)

Visit our Web Site at <http://www.phytotechlab.com>

©2013 PhytoTechnology Laboratories®